


In TrueChem version 21.4.1, clicking the Options button in the Graphs incorrectly positions the Options frame so that it is not accessible via the mouse. In fact, any movement of the mouse immediately closes the Options frame. See figure 1.

This behavior has been corrected starting in version 21.4.2, but if you are using version 21.4.1 you can change a setting in the Options frame by using the Tab key on the keyboard to place the focus to the Options frame. Then use the Up/Down arrow keys on the keyboard to place the focus on the desired setting. Use the space bar to change that item's setting and close the Options frame.

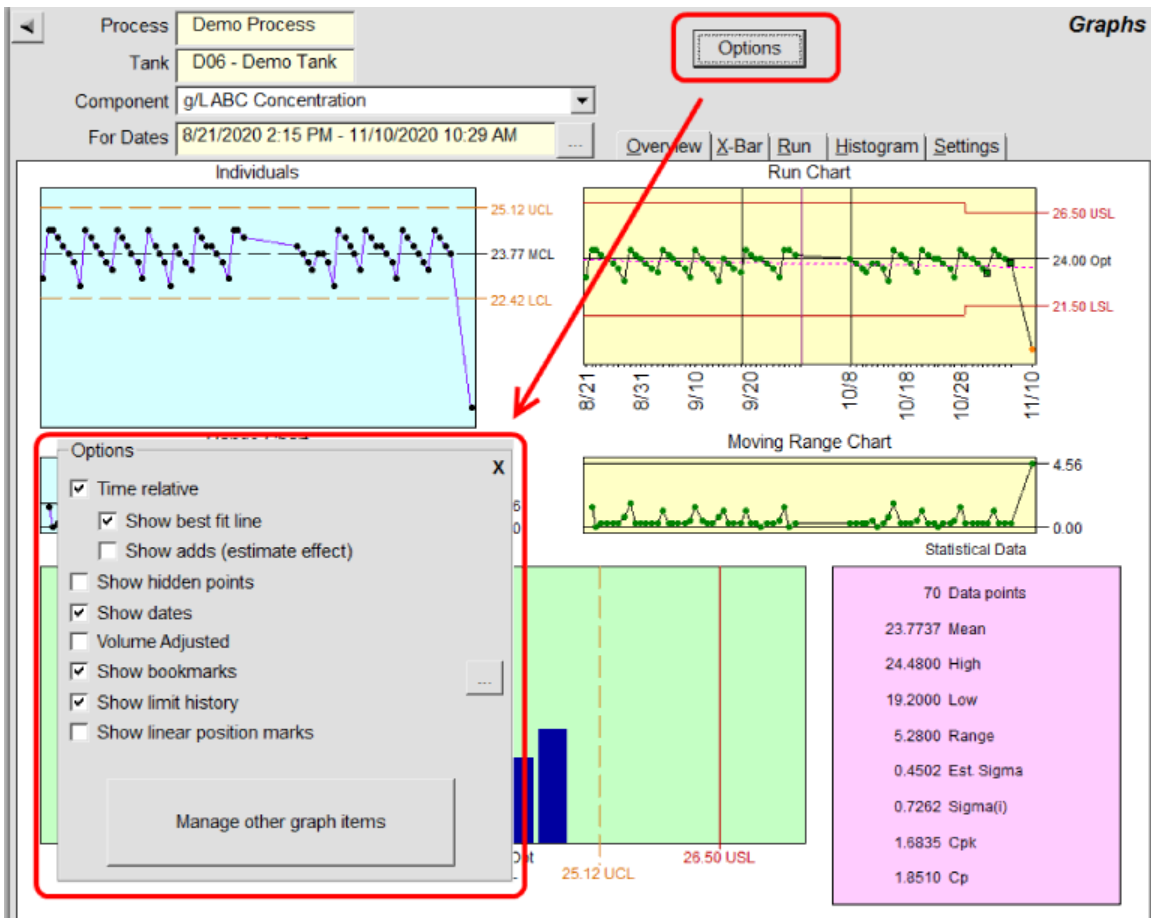

(figure 1)

If you have any questions or need any help, please contact technical support at 1-800-448-2548 ext. 202